

ECOPLUS


LIGHTED 30X MICROSCOPE

With "Rotatable Direct Light Source" Feature

TO OPERATE MICROSCOPE

- 1) Slide out Battery Cover (B), install 2-AA batteries (+) positive end down. Close Battery Cover (Fig. 1).
- 2) Place specimen on flat surface.
- 3) Put microscope over specimen with the Eye Piece (A) directly above specimen.
- 4) Switch on Light Switch (D). For brighter image, turn Bulb Rotation Knob (E) toward side to pinpoint the light beam on specimen.
- 5) Look through eyepiece (A), turn Focusing Wheel (C) to obtain sharpest image.

FIG. 1

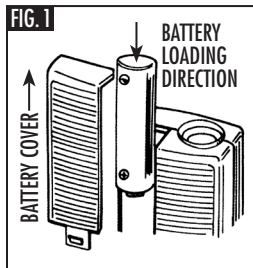

TO REPLACE BULB

- 1) Adjust Bulb (F) upright by turning Bulb Rotation Knob (E) towards center
- 2) Remove Transparent Cover (G) by sliding out (Fig. 2).
- 3) Unscrew faulty bulb (F).
- 4) Replace with new bulb and screw securely.

FIG. 2

